



ZYMO RESEARCH

The Beauty of Science is to Make Things Simple

INSTRUCTION MANUAL

ZR Genomic DNA™-Tissue MiniPrep

Catalog Nos. **D3050 & D3051**

Highlights

- For high quality DNA purification from solid tissues (e.g., tailsnips, earpunches, adipose tissue, etc.), whole blood, plasma, serum, buffy coat, lymphocytes, cultured cells, buccal cells, FFPE tissues, semen, hair, and other biological sources.
- Combines Proteinase K digestion with innovative *Fast-Spin* column purification technology.
- Isolated DNA is ideal for PCR, endonuclease digestion, Southern blotting, bisulfite conversion/methylation detection, sequencing, genotyping, etc.

Contents

Product Contents & Specifications.....	1
Product Description	2
General Considerations.....	3
Reagent Preparation.....	4
Protocols	
Solid Tissue.....	4
Whole Blood, Serum, and Plasma	5
Cell Monolayer.....	5, 6
Biological Liquids/Cell Suspensions	7
Alternative Protocols for Hair, Feathers, and FFPE Tissue....	8
Troubleshooting.....	8
Ordering Information.....	9
Epigenetics-Related Products from Zymo.....	10

Satisfaction of all Zymo Research products is guaranteed. If you are not satisfied with this product please call 1-888-882-9682.

Product Contents

ZR Genomic DNA™-Tissue MiniPrep (Kit Size)	D3050 (50 Preps.)	D3051 (200 Preps.)	Storage Temperature
Proteinase K & Storage Buffer*	2 x 5 mg	2 x 20 mg	-20°C(after mixing)
2X Digestion Buffer**	5 ml	20 ml	Room Temp.
Genomic Lysis Buffer***	50 ml	2 x 100 ml	Room Temp.
DNA Pre-Wash Buffer**	15 ml	50 ml	Room Temp.
g-DNA Wash Buffer	50 ml	100 ml	Room Temp.
DNA Elution Buffer	10 ml	50 ml	Room Temp.
Zymo-Spin™ IIC Columns	50 columns	200 columns	Room Temp.
Collection Tubes	100 tubes	400 tubes	Room Temp.
Instruction Manual	1	1	-

Note - Integrity of kit components is guaranteed for up to one year from date of purchase. Reagents are routinely tested on a lot-to-lot basis to ensure they provide maximal performance and reliability.

* The Proteinase K is stable as shipped. Add 260 µl (1,040 µl for D3051) **Proteinase K Storage Buffer** to each **Proteinase K** tube prior to use. The final concentration of **Proteinase K** after the addition of **Proteinase K Storage Buffer** is ~20 mg/ml.

** The **2X Digestion Buffer** and **DNA Pre-Wash Buffer** may have formed a precipitate. If this is the case, incubate at 37°C to solubilize.

***.For optimal performance, add beta-mercaptoethanol to 0.5%(v/v) i.e., 250 µl per 50 ml or 500 µl per 100 ml.

Specifications

- **Sample Sources** – Solid tissues (e.g., tailsnips, earpunches, adipose tissue, etc.), whole blood, plasma, serum, buffy coat, lymphocytes, cultured cells, buccal cells, FFPE tissues, semen, hair, and other biological sources are effectively processed using this kit.
- **DNA Purity** – High quality DNA for PCR, endonuclease digestion, Southern blotting, bisulfite conversion/methylation detection, sequencing, genotyping, etc., is eluted with **DNA Elution Buffer** or water. ($A_{260}/A_{280} \geq 1.8$)
- **DNA Size** – Capable of recovering genomic and mitochondrial DNA sized fragments from 100 bp to ≥ 40 kb. If present, parasitic, microbial, and viral DNA will also be recovered. Typical fragment sizes range from 25 kb-35 kb.
- **DNA Yield** – The DNA binding capacity of the column is 25 µg. Typically, mammalian tissues yield: 1-3 µg DNA per mg skeletal, heart, and brain tissues and 3-5 µg DNA per mg liver, kidney and lung tissues. Human whole blood will yield 3-7 µg DNA per 100 µl blood sampled. DNA is eluted into ≥ 30 µl **DNA Elution Buffer** or water.
- **Product Detergent Tolerance** – $\leq 5\%$ Triton X-100, $\leq 5\%$ Tween-20, $\leq 5\%$ Sarkosyl, $\leq 0.1\%$ SDS.
- **Equipment** – Water bath or heat block (55°C), microcentrifuge, and vortex.

Note - ™ Trademarks of Zymo Research Corporation. This product is for research use only and should only be used by trained professionals. It is not intended for use in diagnostic procedures. Some reagents included with this kit are irritants. Wear protective gloves and eye protection. Follow the safety guidelines and rules enacted by your research institution or facility.

ZYMO RESEARCH CORP.

Product Description

The **ZR Genomic DNA™-Tissue MiniPrep** is a simple procedure for the rapid isolation of total DNA (e.g., genomic, mitochondrial, parasitic, microbial, viral) from a variety of solid tissues. This product has been optimized for maximal recovery of ultra-pure DNA without RNA contamination and is also compatible with buffy coat, bone marrow, cells from culture, whole blood (fresh or stored), serum, plasma, and many biological liquid samples. For processing, simply digest the sample with the supplied **Proteinase K** then add the **Genomic Lysis Buffer**, vortex, and transfer the mixture to the supplied **Zymo-Spin™ Column**. PCR inhibitors are effectively removed during the purification process and purified DNA is suitable for downstream applications including: PCR, Southern blotting, DNA sequencing, endonuclease digestion, bisulfite conversion/methylation analysis, etc.

The **ZR-96 Genomic DNA™-Tissue MiniPrep** (D3055, D3056, D3057) provides high-throughput (i.e., 96-well plate) processing of solid tissue samples.


Zymo Research offers the following for rapid, precise DNA methylation detection...

1.) **EZ DNA Methylation™ Kit** (D5001, D5002, D5003)

2.) **EZ DNA Methylation-Gold™ Kit** (D5005, D5006, D5007, D5008)


3.) **EZ DNA Methylation-Direct™ Kit** (D5020, D5021, D5022, D5023)

Sample
+
Proteinase K at 55°C

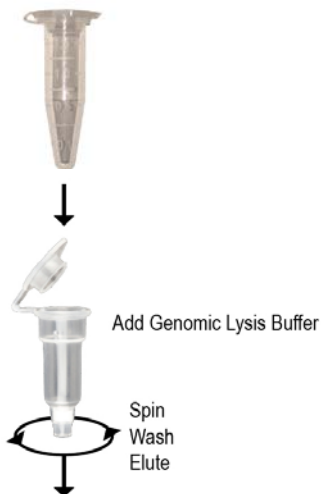

Ultra-pure DNA for...

- ✓ PCR
- ✓ Endonuclease Digestion
- ✓ Southern Blotting
- ✓ Genotyping
- ✓ Bisulfite Conversion & Methylation Analysis

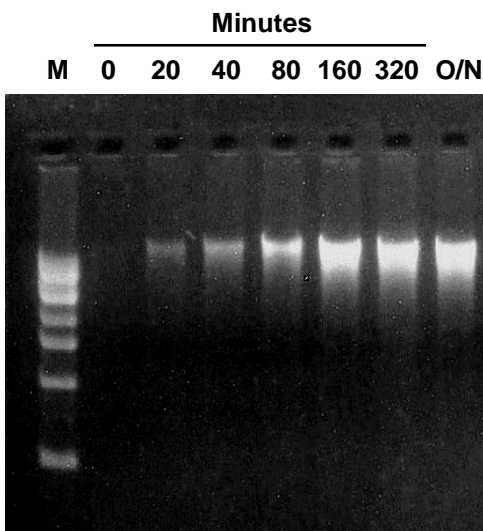

High yield/quality DNA is successfully isolated from porcine muscle using the **ZR Genomic DNA™-Tissue MiniPrep**. Equivalent amounts (25 mg) of muscle tissue were processed using the **ZR Genomic DNA™-Tissue MiniPrep** after incubation with Proteinase K at 55°C for the indicated times (in minutes) or overnight (O/N). Equal volumes of eluted DNA were then analyzed in a 0.8% (w/v) TAE/agarose/ethidium bromide gel. The size marker "M" is a 1 kb ladder (Zymo Research).

Please visit:
www.zymoresearch.com
 for a comprehensive list of
 genomic DNA purification
 products.

General Considerations When Purifying Genomic DNA

Zymo Research offers a range of genomic DNA isolation kits that are suitable for extracting high molecular weight DNA from a wide variety of sample types. Kits are tailor-made for specific applications and feature chemical, Proteinase K, and/or mechanical lysis technologies depending on the starting material (see table below).

DNA Extraction Method	Applications
Chemical	<i>Soft tissue samples</i> from humans, mice, etc., including: whole blood, plasma, serum, cells, buffy coat, buccal cells, biological liquids, crude homogenates, etc.
Proteinase K & Chemical	<i>Solid tissue samples</i> from humans, mice, etc., including: tailsnips, earpunches, hair*, feathers*, and FFPE* samples, as well as all of the above.
Mechanical Homogenization & Chemical	<i>Tough tissues and organisms</i> including: insects, arthropods, fungi, gram (+/-) bacteria, and microorganisms in soil, sludge, feces, or water, as well as most of the above.

The **ZR Genomic DNA™-Tissue MiniPrep** includes Proteinase K digestion and chemical lysis for the rapid, efficient purification of DNA (up to 25 µg/prep.) from soft and solid tissues, cells, and a range of biological liquids (see table below for sample types and protocol recommendations).

Recommended Protocol	Sample Types
Solid Tissue	<i>Solid tissue samples</i> from humans, mice, etc., including: tailsnips, earpunches, hair*, feathers*, and FFPE* samples. (pg. 4)
Whole Blood, Serum, and Plasma	Whole blood, plasma, and serum. (pg. 5)
Cell Monolayer	Monolayer cells ($\leq 5 \times 10^6$) from culture. (pgs. 5-6)
Biological Liquids and Cell Suspensions	Biological liquids including: semen, CSF, buffy coat, body fluids. Cell suspensions containing less than 5×10^6 cells (e.g., buffy coat, suspension cultured cells, etc.) (pg. 7)

* With protocol modification. See Alternative Protocols **(pg. 8.)**

Starting Material: The quality of the sampled material will affect both the yield and quality of the purified DNA. Freshly sampled tissues and cells yield the highest quantity/ quality DNA. If sampling from “stored” sources and/or if samples have been subject to repeated freeze/thawing, yields may decrease and the purified DNA may be degraded (e.g., FFPE).

Removal of PCR Inhibitors: The **ZR Genomic DNA™-Tissue MiniPrep** has been designed for the efficient removal of PCR inhibitors during DNA purification from the samples listed in the tables above. However, some environmental samples including soil, plants, and manure (feces) will require alternative technologies (see sidebar) for the effective removal of polyphenolic PCR inhibitors.

Storage of Purified DNA: The eluted DNA can be used immediately for molecular-based applications or stored $\leq -20^\circ\text{C}$.

The **ZR Soil Microbe DNA Kit™** (D6001), **ZR Fecal DNA Kit™** (D6010), and **ZR Plant/Seed DNA Kit™** (D6020) can be used for the purification of inhibitor-free DNA from soil, feces, and plants, respectively.

ZYMO RESEARCH CORP.

Reagent Preparation

- Add 260 μl (1,040 μl for D3051) **Proteinase K Storage Buffer** to each **Proteinase K** tube prior to use. The final concentration of **Proteinase K** after the addition of **Proteinase K Storage Buffer** is ~20 mg/ml.
- For optimal performance, add beta-mercaptoethanol (user supplied) to the **Genomic Lysis Buffer** to a final dilution of 0.5%(v/v) i.e., 250 μl per 50 ml or 500 μl per 100 ml.

Protocols

Solid Tissue – Including: tailsnips, earpunches, biopsies, etc.

The following is for the purification of DNA from up to 25 mg fresh or frozen tissue. Typical yields are: 1-3 μg DNA per mg skeletal, heart, and brain tissues and 3-5 μg DNA per mg liver, kidney, and lung tissues. For hair, feathers and FFPE tissues follow **Alternative Protocols I** and **II** on page 8, respectively.

1. To a tissue sample (≤ 25 mg) in a microcentrifuge tube add a solution of...

H ₂ O	95 μl
2X Digestion Buffer	95 μl
Proteinase K	10 μl

2. Mix and then incubate the tube at 55°C for 1-3 hours.

Note: If required (e.g., FFPE samples), digesting samples overnight at 55°C with Proteinase K is possible without affecting the integrity of the DNA.

3. Add 700 μl **Genomic Lysis Buffer** to the tube and mix thoroughly by vortexing. Centrifuge at 10,000 x *g* for one minute to remove insoluble debris.
4. Transfer the supernatant to a **Zymo-Spin™ IIC Column** in a **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
5. Add 200 μl of **DNA Pre-Wash Buffer** to the spin column in a new **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
6. Add 400 μl of **g-DNA Wash Buffer** to the spin column. Centrifuge at 10,000 x *g* for one minute.
7. Transfer the spin column to a clean microcentrifuge tube. Add ≥ 50 μl **DNA Elution Buffer** or water (e.g., add 200 μl if sampling 25 mg tissue) to the spin column. Incubate 2-5 minutes at room temperature, then centrifuge at top speed for 30 seconds to elute the DNA. The eluted DNA can be used immediately for molecular based applications or stored $\leq -20^\circ\text{C}$ for future use.

For **Technical Assistance**, please contact 1-888-882-9682 or E-mail tech@zymoresearch.com.

Incubate 12-16 hours for formalin-fixed deparaffinized samples.

The column capacity is ~1 ml.

Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0 . Also, the total yield may be improved by eluting the DNA with Elution Buffer or water pre-equilibrated to 60-70°C or by performing and pooling sequential elutions.

Human whole blood should yield between 3-7 µg DNA per 100 µl.

The column capacity is ~1 ml.

Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0. Also, the total yield may be improved by eluting the DNA with Elution Buffer or water pre-equilibrated to 60-70°C or by performing and pooling sequential elutions.

Generally, no more than 5×10^6 cells should be sampled, for larger samples will exceed the binding capacity of the spin column.

Whole Blood, Serum and Plasma

The following is for the purification of DNA from up to 100 µl whole blood, serum or plasma (the volumes can be adjusted depending on your requirements). Fresh, frozen, or preserved blood (in EDTA, citrate, or heparin) can be used.

1. Adjust total volume of sample (blood, serum, or plasma) to 100 µl with water in a microcentrifuge tube and then add the following...

2X Digestion Buffer	95 µl
Proteinase K	5 µl

Example: Add 40 µl H₂O to 60 µl blood, serum, or plasma prior to adding the 2X Digestion Buffer and Proteinase K.

2. Mix and then incubate the tube at 55°C for 20 minutes.
3. Add 700 µl **Genomic Lysis Buffer** to the tube and mix thoroughly by vortexing.
4. Transfer the mixture to a **Zymo-Spin™ IIC Column** in a **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
5. Add 200 µl of **DNA Pre-Wash Buffer** to the spin column in a new **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
6. Add 400 µl of **g-DNA Wash Buffer** to the spin column. Centrifuge at 10,000 x *g* for one minute.
7. Transfer the spin column to a clean microcentrifuge tube. Add ≥50 µl **DNA Elution Buffer** or water (e.g., add 100 µl if sampling 100 µl blood, serum, or plasma) to the spin column. Incubate 2-5 minutes at room temperature, then centrifuge at top speed for 30 seconds to elute the DNA. The eluted DNA can be used immediately for molecular based applications or stored ≤-20°C for future use.

Cell Monolayer

The following procedure is designed for up to 5×10^6 monolayer cells. Although cell types and culture conditions may vary, the protocol will work with high-density growth cells (e.g., HeLa cells) as well as with low-density growth cells (e.g., neuronal cells).

1. Trypsinize or scrape adherent cells from a culture flask or plate. Centrifuge the suspension at approximately 500 x *g* for 5 minutes. Remove the supernatant and resuspend the cell pellet in 1 ml PBS (Phosphate Buffered Saline) and then transfer suspension to a microcentrifuge tube. Centrifuge the suspension at approximately 500 x *g* for 5 minutes. Remove the supernatant and resuspend the pellet in a solution of...

H ₂ O	95 µl
2X Digestion Buffer	95 µl
Proteinase K	5 µl

(continued on next page)

ZYMO RESEARCH CORP.

2. Incubate the tube at 55°C for 20 minutes.
3. Add 700 µl **Genomic Lysis Buffer** to the tube and mix thoroughly by vortexing. Centrifuge at 10,000 x *g* for one minute to remove insoluble debris.
4. Transfer the supernatant to a **Zymo-Spin™ IIC Column** in a **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
5. Add 200 µl of **DNA Pre-Wash Buffer** to the spin column in a new **Collection Tube**. Centrifuge at 10,000 x *g* for one minute.
6. Add 400 µl of **g-DNA Wash Buffer** to the spin column. Centrifuge at 10,000 x *g* for one minute.
7. Transfer the spin column to a clean microcentrifuge tube. Add ≥50 µl **DNA Elution Buffer** or water (e.g., add 200 µl if sampling 5x10⁶ cells) to the spin column. Incubate 2-5 minutes at room temperature, then centrifuge at top speed for 30 seconds to elute the DNA. The eluted DNA can be used immediately for molecular based applications or stored ≤-20°C for future use.

The column capacity is ~1 ml.

Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0. Also, the total yield may be improved by eluting the DNA with Elution Buffer or water pre-equilibrated to 60-70°C or by performing and pooling sequential elutions.

Guidelines for Monolayer Cell DNA Isolation: Cell numbers (growth densities) can vary between different cell types. Table 1 (below) provides an approximation of the cell numbers that can be recovered from different culture containers for “high-density” growth cells like CV1 and HeLa cells.

Table 1: Culture Plate/Flask Growth Area (cm²) and Cell Number

Culture Container	Well /Flask Surface Area	Cell Number
96-well plate	0.32-0.6 cm ²	4-5x10 ⁴
24-well plate	2 cm ²	1-3x10 ⁵
12-well plate	4 cm ²	4-5x10 ⁵
6-well plate	9.5 cm ²	0.5-1x10 ⁶
T25 Culture Flask	25 cm ²	2-3x10 ⁶
T75 Culture Flask	75 cm ²	0.6-1x10 ⁷
T175 Culture Flask	175 cm ²	2-3x10 ⁷

Biological Liquids and Cell Suspensions

Cells should be processed directly from biological fluids or from suspension in PBS, TE, or compatible buffers.

The following protocol is designed for up to 100 µl of biological liquid sample including semen, CSF, buffy coat, body fluids, and cell suspensions containing less than 5×10^6 cells.

1. Adjust total volume of liquid sample to 100 µl with water in a microcentrifuge tube and then add the following...

2X Digestion Buffer	95 µl
Proteinase K	5 µl

2. Mix and then incubate the tube at 55°C for 20 minutes.
3. Add 700 µl **Genomic Lysis Buffer** to the tube and mix thoroughly by vortexing.
4. Transfer the mixture to a **Zymo-Spin™ IIC Column** in a **Collection Tube**. Centrifuge at 10,000 x g for one minute.
5. Add 200 µl of **DNA Pre-Wash Buffer** to the spin column in a new **Collection Tube**. Centrifuge at 10,000 x g for one minute.
6. Add 400 µl of **g-DNA Wash Buffer** to the spin column. Centrifuge at 10,000 x g for one minute.
7. Transfer the spin column to a clean microcentrifuge tube. Add ≥50 µl **DNA Elution Buffer** or water (e.g., add 200 µl if sampling liquids containing 5×10^6 cells) to the spin column. Incubate 2-5 minutes at room temperature, then centrifuge at top speed for 30 seconds to elute the DNA. The eluted DNA can be used immediately for molecular based applications or stored ≤-20°C for future use.

The column capacity is ~1 ml.

Elution of DNA from the column is dependent on pH and temperature. If water is used, ensure the pH is >6.0. Also, the total yield may be improved by eluting the DNA with Elution Buffer or water pre-equilibrated to 60-70°C or by performing and pooling sequential elutions.

Alternative Protocols:

I.) For Hair, Feathers, or Related Samples: Freshly prepared DTT (dithiothreitol) (not provided) needs to be added to Step 1 of the Solid Tissue Protocol (page 3) as follows...

H ₂ O	90 µl
2x Digestion Buffer	90 µl
DTT (1 M)	10 µl
Proteinase K	10 µl

Then follow with the rest of the procedure as indicated.

II.) For FFPE Samples: Tissues need to be deparaffinized prior to Step 1 of the Solid Tissue Protocol (page 4) by...

- i. Removing (trimming) as much paraffin from the sample(s) as possible.
- ii. Transfer samples to 1.5 ml microcentrifuge tubes. Add 750 µl xylene (not provided) to the samples.
- iii. Vortex and incubate samples at room temperature for 1 hour with gentle rocking.
- iv. Centrifuge for 1 minute at 10,000 x g and remove the xylene from the sample. Repeat steps 2-4.
- v. Wash two times with 1 ml EtOH (100%) for 5 minutes with gentle rocking.
- vi. Wash two times with 1 ml EtOH (95%) for 5 minutes with gentle rocking.
- vii. Wash two times with 1 ml EtOH (75%) for 5 minutes with gentle rocking.
- viii. Wash once with 1 ml ddH₂O for 5 minutes with gentle rocking. Remove as much water from the sample as possible
- ix. Use sample or store at -80°C.

Note: For steps v-viii, add the wash, vortex briefly, and incubate for 5 minutes with gentle rocking. Remove wash from the sample after centrifugation at 10,000 x g for 1 minute.

Troubleshooting:

1. DNA degradation: Check for DNase contamination. All reagents and components supplied with the **ZR Genomic DNA™-Tissue MiniPrep** are DNase-free. However, DNase contamination can result during the processing of some samples. Check pipets, pipet tips, microcentrifuge tubes, etc., and exercise the appropriate precautions during the DNA purification procedure. Make sure Proteinase K digestions are performed at 55°C as indicated.
2. DNA is not performing well in subsequent experiments: Ensure the correct volume of **Genomic Lysis Buffer** has been added to the sample. Also, make sure all centrifugation steps are completed for the indicated times and speeds (rcfs). Failure to do so may result in incomplete washing, which may cause salts to be eluted with the DNA affecting quantitation and subsequent experiments including enzymatic processes like PCR.
3. RNA contamination: The buffers and spin columns provided in this kit are designed to efficiently remove RNA during the DNA purification procedure. However, additional RNA removal (e.g., digestion with RNase A) may be necessary for subsequent applications sensitive to trace amounts of RNA.

Ordering Information

Product Description	Catalog No.	Kit Size
ZR Genomic DNA™-Tissue MicroPrep	D3040	50 preps.
	D3041	200 preps.
ZR Genomic DNA™-Tissue MiniPrep	D3050	50 preps.
	D3051	200 preps.
ZR Genomic DNA™-Tissue MidiPrep	D3110	25 preps.
ZR-96 Genomic DNA™-Tissue MiniPrep	D3055	2x96 preps.
	D3056	4x96 preps.
	D3057	10x96 preps.

For Individual Sale	Catalog No.	Amount
Proteinase K & Storage Buffer	D3001-2-5	5 mg set
	D3001-2-20	20 mg set
2X Digestion Buffer	D3050-1-5	5 ml
	D3050-1-20	20 ml
Genomic Lysis Buffer	D3004-1-50	50 ml
	D3004-1-100	100 ml
DNA Pre-Wash Buffer	D3004-5-15	15 ml
	D3004-5-30	30 ml
	D3004-5-50	50 ml
g-DNA Wash Buffer	D3004-2-50	50 ml
	D3004-2-100	100 ml
DNA Elution Buffer	D3004-4-4	4 ml
	D3004-4-10	10 ml
	D3004-4-50	50 ml
Zymo-Spin™ IIC Columns	C1011-50	50 columns
	C1011-250	250 columns
Collection Tubes	C1001-50	50 tubes
	C1001-500	500 tubes
	C1001-1000	1,000 tubes

ZYMO RESEARCH CORP.

Popular DNA Purification Products from Zymo Research

Product	Format	Kit Size	Cat No.
Fragment DNA Clean-up, Concentration & Recovery			
DNA Clean & Concentrator™-5	Spin Column Format (up to 5 µg/prep.)	50 preps. 200 preps.	D4003*, D4013 D4004*, D4014
DNA Clean & Concentrator™-25	Spin Column Format (up to 25 µg/prep.)	50 preps. 200 preps.	D4005*, D4033 D4006*, D4034
DNA Clean & Concentrator™-100	Spin Column Format (up to 100 µg/prep.)	25 preps. 50 preps.	D4029 D4030
DNA Clean & Concentrator™-500	Spin Column Format (up to 500 µg/prep.)	10 preps. 20 preps.	D4031 D4032
ZR-96 DNA Clean & Concentrator™-5	96-Well Format (up to 5 µg/well; deep well)	2x96 preps. 4x96 preps.	D4023 D4024
Genomic DNA Clean & Concentrator™	Spin Column Format (up to 10 µg/prep.)	25 preps. 100 preps.	D4010 D4011
ZR-96 DNA Clean-up Kit™	96-Well Format (up to 5 µg/well; shallow well)	2x96 preps. 4x96 preps.	D4017 D4018
ZR DNA Sequencing Clean-up Kit™	Spin Column Format (up to 5 µg/prep.)	50 preps. 200 preps.	D4050 D4051
ZR-96 DNA Sequencing Clean-up Kit™	96-Well Format (up to 5 µg/well)	2x96 preps. 4x96 preps.	D4052 D4053
OneStep™ PCR Inhibitor Removal Kit	Spin Column Format (up to 25 µg/prep.)	50 preps.	D6030
OneStep-96™ PCR Inhibitor Removal Kit	96-Well Format (up to 5 µg/well)	2x96 preps.	D6035
Zymoclean™ Gel DNA Recovery Kit	Spin Column Format (up to 5 µg/prep.)	50 preps. 200 preps.	D4001 D4002
ZR-96 Zymoclean™ Gel DNA Recovery Kit	96-Well Format (up to 5 µg/well)	2x96 preps. 4x96 preps.	D4021 D4022
Zymoclean™ Large Fragment DNA Recovery Kit	Spin Column Format (up to 10 µg/prep.)	25 preps. 100 preps.	D4045 D4046
Plasmid DNA Isolation			
Zyppy™ Plasmid Miniprep Kit	Pellet Free, Spin Column Format	50 preps. 100 preps. 400 preps. 800 preps.	D4036 D4019 D4020 D4037
Zyppy™ Plasmid Midiprep Kit	Pellet Free, Spin Column Format	25 preps. 50 preps.	D4025 D4026
Zyppy™ Plasmid Maxiprep Kit	Spin/Vacuum Column Format	10 preps. 20 preps.	D4027 D4028
ZR Plasmid Miniprep™-Classic	Spin Column Format	100 preps. 400 preps. 800 preps.	D4015 D4016 D4054
ZR BAC DNA Miniprep Kit	BAC/PAC plasmid DNA Isolation. Spin Column Format	25 preps. 100 preps.	D4048 D4049
Environmental DNA Isolation			
ZR Soil Microbe DNA MicroPrep™	Bead Bashing, Spin Column Format (up to 5 µg/prep.)	50 preps.	D6003
ZR Soil Microbe DNA MiniPrep™	Bead Bashing, Spin Column Format (up to 25 µg/prep.)	50 preps.	D6001
ZR Soil Microbe DNA MidiPrep™	Bead Bashing, Spin Column Format (up to 125 µg/prep.)	25 preps.	D6101
ZR-96 Soil Microbe DNA Kit™	Bead Bashing, 96-Well Format (up to 5 µg/well)	2x96 preps.	D6002
ZR Fungal/Bacterial DNA MicroPrep™	Bead Bashing, Spin Column Format (up to 5 µg/prep.)	50 preps.	D6007
ZR Fungal/Bacterial DNA MiniPrep™	Bead Bashing, Spin Column Format (up to 25 µg/prep.)	50 preps.	D6005
ZR Fungal/Bacterial DNA MidiPrep™	Bead Bashing, Spin Column Format (up to 125 µg/prep.)	25 preps.	D6105
ZR-96 Fungal/Bacterial DNA Kit™	Bead Bashing, 96-Well Format (up to 5 µg/well)	2x96 preps.	D6006
ZR Fecal DNA MicroPrep™	Bead Bashing, Spin Column Format (up to 5 µg/prep.)	50 preps.	D6012
ZR Fecal DNA MiniPrep™	Bead Bashing, Spin Column Format (up to 25 µg/prep.)	50 preps.	D6010
ZR Fecal DNA MidiPrep™	Bead Bashing, Spin Column Format (up to 125 µg/prep.)	25 preps.	D6110
ZR-96 Fecal DNA Kit™	Bead Bashing, 96-Well Format (up to 5 µg/well)	2x96 preps.	D6011
ZR Tissue & Insect DNA MicroPrep™	Bead Bashing, Spin Column Format (up to 5 µg/prep.)	50 preps.	D6015
ZR Tissue & Insect DNA MiniPrep™	Bead Bashing, Spin Column Format (up to 25 µg/prep.)	50 preps.	D6016
ZR Tissue & Insect DNA MidiPrep™	Bead Bashing, Spin Column Format (up to 125 µg/prep.)	25 preps.	D6115
ZR-96 Tissue & Insect DNA Kit™	Bead Bashing, 96-Well Format (up to 5 µg/well)	2x96 preps.	D6017
ZR Plant/Seed DNA MicroPrep™	Bead Bashing, Spin Column Format (up to 5 µg/prep.)	50 preps.	D6022
ZR Plant/Seed DNA MiniPrep™	Bead Bashing, Spin Column Format (up to 25 µg/prep.)	50 preps.	D6020
ZR Plant/Seed DNA MidiPrep™	Bead Bashing, Spin Column Format (up to 125 µg/prep.)	25 preps.	D6120
ZR-96 Plant/Seed DNA Kit™	Bead Bashing, 96-Well Format (up to 5 µg/well)	2x96 preps.	D6021

* Uncapped Spin Column Format

ZYMO RESEARCH CORP.

Phone: (949) 679-1190 ▪ Toll Free: (888) 882-9682 ▪ Fax: (949) 266-9452 ▪ info@zymoresearch.com ▪ www.zymoresearch.com